

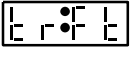
# Foot Switch

## Part #: ADPT-FOOTSW-2

### Description

The Foot Switch cable allows the operator to start and stop a syringe pump remotely by pressing the switch with their foot.

### Pump Setup

- Set the “Operational Trigger”, of the pump to “foot switch”. To set, with the pump stopped, press and hold the “Diameter” key to access the pump setup. Wait until the display shows the current “Operational Trigger” [tr:aa]. Use the right-most arrow key to select the mode.  
“foot switch (Ft)” mode is selected when the display shows: . The settings will be stored in non-volatile memory after the display blinks.
- Before connecting the cable, turn power off to the pump.
- Connect the Foot Switch cable connector to the connector on the back of the pump labeled “TTL-I/O”. Secure the cable to the pump with the 2 screws on the cable connector.
- When the power is turned on to the pump, it is ready for use. While the Foot Switch is attached, the 'Start/Stop' button on the front panel can still be used.

