

RS-232 Pump Network Secondary Cable

<u>Part #</u>	<u>Description</u>
CBL-NET-7	7' Pump network secondary cable
CBL-NET-25	25' Pump network secondary cable

Contents:

- 1) Pump to pump network cable, 7' or 25' (Figure 1).



Figure 1: Network Cable (7' or 25')

Instructions:

- 1) Attach either end of the "Network cable" to the **"To Network"** RJ-11 (telephone style) connector on the back of the previous pump on the pump network.
- 2) Attach the other end of the "Network cable" to the **"To Computer"** RJ-11 (telephone style) connector of the next pump on the pump network
- 3) Refer to the pump's User's Manual to setup the baud rate and network address. The computer and each pump must have the same baud rate.
 - Each pump on the pump network must have a unique network address.

NOTE: The first pump is attached to the computer using cable CBL-PC-PUMP.

